



Zettachring 6, 70567 Stuttgart, Germany  
Tel: +49-711-13257-0 Fax: +49-711-13257-89  
E-Mail: info.sep.eu@tosoh.com  
Web: www.tosohbioscience.com

3604 Horizon Drive, Suite 100, King of Prussia, PA 19406, USA  
Tel: +1 800-366-4875 Fax: +1 610-272-3028  
E-Mail: info.sep.am@tosoh.com  
Web: www.tosohbioscience.com

## OPERATING CONDITIONS and SPECIFICATIONS

### TSK-GEL® G6000PW Products

<b>Part Numbers:</b>	05765	7.5 mm ID x 30.0 cm L	17 µm
	05109	7.5 mm ID x 60.0 cm L	17 µm
	05154	21.5 mm ID x 60.0 cm L	23 µm

This sheet contains the recommended operating conditions and the specifications for TSK-GEL G6000PW columns. Installation instructions and column care information are described in a separate Instruction Manual.

#### A. OPERATING CONDITIONS

- Shipping Solvent: Water
- Max. Flow Rate: 1.2 ml/min 7.5mm ID  
8.0 ml/min 21.5mm ID

**NOTE:** When a buffer with high viscosity is used, the maximum flow rate may have to be reduced so as not to exceed the maximum pressure drop. When changing solvents, use a flow rate equal to 25% of the maximum flow rate.

- Standard Flow Rate: 0.5 - 1.0 ml/min 7.5mm ID  
1.6 - 6.0 ml/min 21.5mm ID
- Max. Pressure: 10 kg/cm<sup>2</sup> = 150 psi 7.5mm ID x 30.0 cm L and 21.5mm ID x 60.0 cm L  
20 kg/cm<sup>2</sup> = 300 psi 7.5mm ID x 60.0 cm L
- pH Range: 2.0 - 12.0
- Salt Conc.: ≤ 0.5 Molar
- Organic Conc.: ≤ 20%
- Temperature: 10 - 80°C Reduce flow rate when operating below 10°C.

- Cleaning Solvents:
  - (1) High salt concentration buffer (0.5 - 1.0 M), or
  - (2) pH 2 - 3 or pH 9 - 12 buffer, or
  - (3) Buffer with acetonitrile or methanol, or if nothing else is successful
  - (4) Buffer with urea or SDS

**NOTE:** Choose a cleaning solvent based on sample properties, e.g. use (1) to remove basic polymers, and (3) to remove hydrophobic proteins etc.

- Storage: Store the column in a 0.05% NaN<sub>3</sub> solution or 20% ethanol in DI water when it will not be used the next day. For overnight storage flush the column at low flow rate with the mobile phase. Prevent air from entering the column!
- Column Protection: The use of a guard column (TSKgel PW Guard Column P/N 06762 for 7.5 m ID or P/N 06758 for 21.5 mm ID) is recommended to prolong the life of the analytical column. Guard column life depends greatly on sample cleanliness. As a general rule, guard columns should be replaced after every 30-40 sample injections, when the peaks become excessively wide, or when the peaks show splitting.

#### B. SPECIFICATIONS

The performance of TSK-GEL G6000PW columns is tested under the conditions described in the Data Sheet. All columns have passed the following quality control specifications

Number of Theoretical Plates (N): ≥ 3,000 7.5mm ID x 30.0 cm L  
≥ 6,000 60.0 cm L columns

Asymmetry Factor (AF): 0.7 - 1.6